



17.02 / -932L+ → -932L+ / 59.00
 17.02 / -932L- → -932L- / 59.00
 43.04 / -PE103 → -PE103 / 74.00

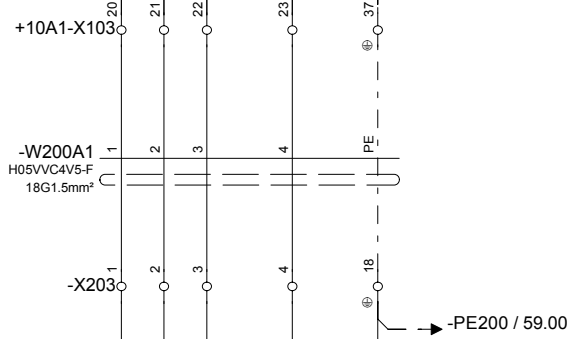


-U2001
53.00
WAGO
Profibus Interface
750-301

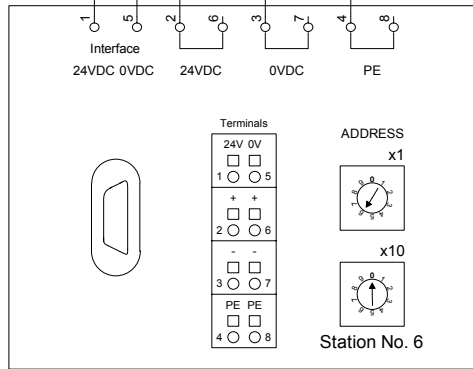

-20A1 []

=FD+20A1/53										=FD+20A1/55	
A	Subsupplier	21.08.2003	G. Grob	Date	03.04.2003	Novo Nordisk A/S 2820 Gentofte/Denmark Niels Steensvej 1/Building HAA	STERIS STERIS GmbH Kalscheuren Str. 02 50354 Hürth Deutschland	-U2001 Profibus DP interface	Freeze-Dryer	FCM 30-D/LN2	= FD + 20A1
Rev.	Modification	Date	Name	Norm	Origin	Crea. f.	Crea. by	Copyright reserved	1383-01	Drawing-No. 1383-01-1001	Page 54 184 Pg

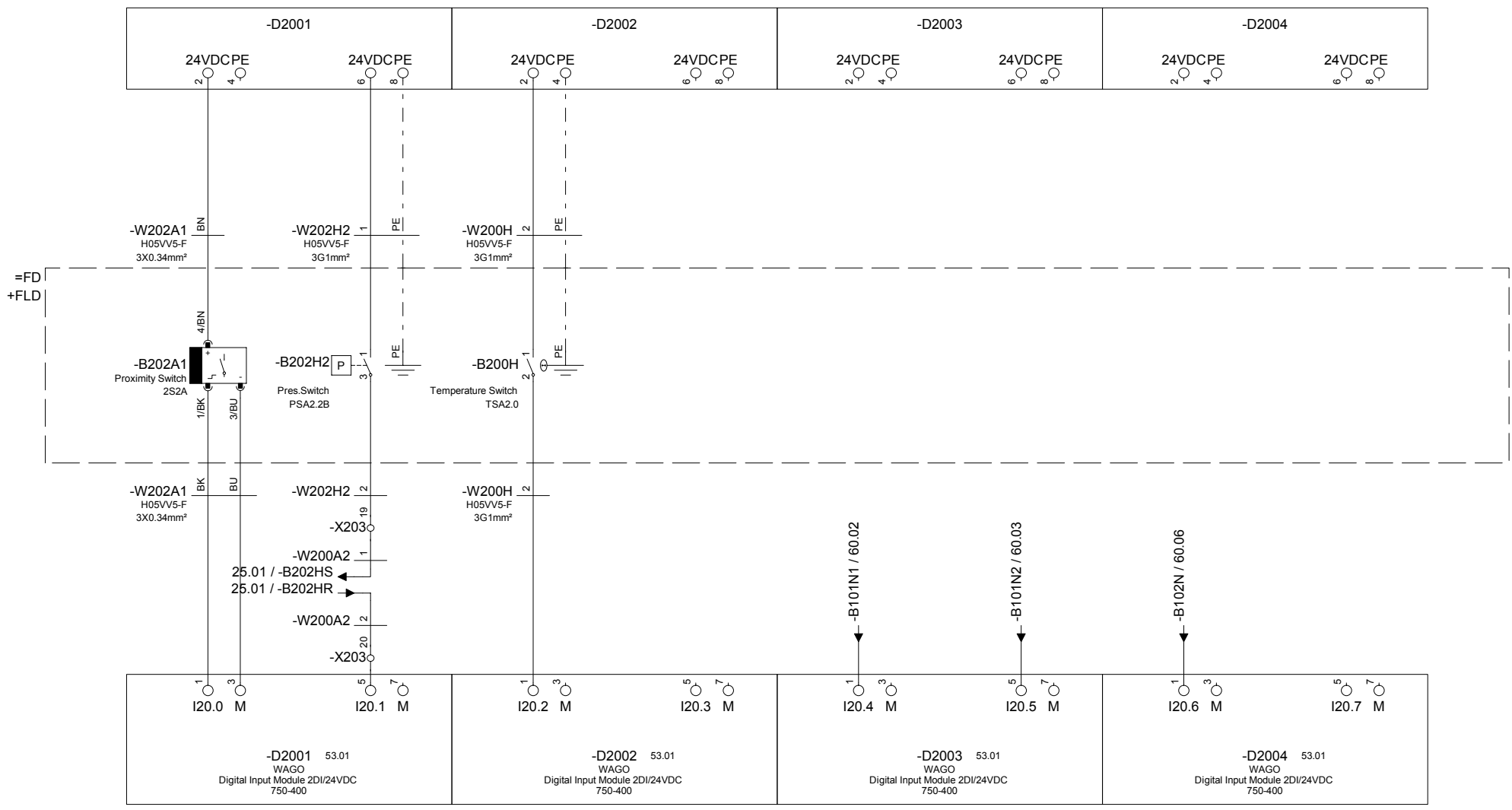

- | | | | | | | | |
|--|---|---|-------|---|---|--|-------|
| Proximity Switch
-B202A1 (2S2A)
Chamber/Condenser Valve
-Y202B (2V2A)
open | Pressure Switch
-B202H2 (PSA 2.2B)
Pressure at Valve
-Y202R (2V2B)
Above 57 Bar | Temperature Switch
-B200H (TSA 2.0)
Hydraulic Oil
Excess Temperature | Spare | Pressure Transmitter
-B101N1 (PIRCA 1.1A)
Chamber Pressure
Vacuum
Heating Element
No Malfunction | Pressure Transmitter
-B101N2 (PIRCA 1.1B)
Chamber Pressure
Vacuum
Heating Element
No Malfunction | Pressure Transmitter
-B102N (PIRA 1.2)
Condenser Pressure
Vacuum
Heating Element
No Malfunction | Spare |
|--|---|---|-------|---|---|--|-------|